



Optical Cryostat - Large Sample Space

The CS202I-DMX-4SS provides the same high performance as our CS202I-DMX-1SS with its all welded stainless steel construction and welded stainless steel instrumentation skirt but provides a larger than standard sample space. The vacuum shroud comes standard with 4 window ports, however a 5th port can be added on the end. The system is capable of vacuum levels of 10^{-7} Torr with an appropriate vacuum pump.

Applications

- Large Samples
- Optical
- Raman
- UV, VIS, IR
- FTIR
- Electro & Photoluminescence
- Resistivity/Hall Probe Experiments
- Diamond Anvil Cell
- Magneto-Optical
- PITS / DLTS
- Thermal, Electrical and Magnetic Susceptibility
- Magneto Optical Kerr Effect (MOKE)

Features

- Cryogen Free, Low Power
- High Performance Stainless Steel Construction
- Large clear view optical windows (1.5 in)
- Large sample viewing angle for optical collection (F/1.1)
- Can operate in any orientation
- Fully customizable

Typical Configuration

- Cold head (DE-202Al)
- Compressor (ARS-4HW)
- 2 Helium Hoses
- Stainless Steel vacuum shroud with 4 window ports for optical and electrical measures (DMX-4SS)
- Nickel Plated OFHC radiation shield
- 2 High purity quartz windows
- Instrumentation for temperature measurement and control:
 - 10 pin hermetic feed through
 - 36 ohm thermfoil heater
 - Silicon diode sensor curve matched to ($\pm 0.5K$) for control
 - Calibrated silicon diode sensor (± 12 mk) with 4 in. free length for accurate sample measurement.
- Wiring for electrical experiments:
 - 10 pin hermetic feed through
 - 4 copper wires
- Sample holder for optical and electrical experiments
- Temperature Controller

Options and Upgrades

- 4K Coldhead (0.1W @ 4.2K)
- 5.5K Coldhead (1W @ 10K)
- 450K High Temperature Interface
- 800K High Temperature Interface
- Turbo upgrade for faster cooldown times
- Custom temperature sensor configuration (please contact our sales staff)
- Custom wiring configurations (please contact our sales staff)
- Window material upgrades (custom materials available)
- Sample holder upgrades (custom sample holders available)



The above picture shows a cryocooler with a vacuum shroud, radiation shield, and sample holder installed.



The above picture shows a complete system (minus the vacuum pump and temperature controller).



Optical Cryostat - Large Sample Space

Cooling Technology

DE-202	Closed Cycle Cryocooler
Refrigeration Type	Pneumatically Driven GM Cycle
Liquid Cryogen Usage	None, Cryogen Free

Temperature*

DE-202AI	< 10K - 350K
DE-202PI	< 5.5K - 350K
DE-202SI	< 4.2K - 350K
With 800K Interface	(Base Temp + 2K) - 700K
With 450K Interface	(Base Temp + 2K) - 450K
Stability	0.1K
*Based on bare cold head with a closed radiation shield, and no additional sources of experimental or parasitic heat load	

Sample Space

Diameter	47.5 mm (1.87in.)
Height	90 mm (3.54 in.)
Sample Holder Attachment	1/4 - 28 screw
Sample Holder	www.arscryo.com/Products/SampleHolders.html

Optical Access

Window Ports	4 - 90° Apart
Diameter	51 mm (2 in)
Clear View	38 mm (1.5 in)
#/F	1.1
Window Material	www.arscryo.com/Products/WindowMaterials.html

Temperature Instrumentation and Control (Standard)

Heater	36 ohm Thermofoil Heater anchored to the coldtip
Control Sensor	Curve Matched Silicon Diode installed on the coldtip
Sample Sensor	Calibrated Silicon Diode with free length wires

Contact ARS for other options

Instrumentation Access

Instrumentation Skirt	Welded Stainless Steel
Pump out Port	1 - NW 25
Instrumentation Ports	3
Instrumentation Wiring	Contact sales staff for options

Vacuum Shroud

Material	Stainless Steel
Length	378 mm (14.9 in)
Diameter	95mm (3.75 in) (at the sample space)
Width	95 mm (3.75 in) (at the sample space)

Radiation Shield

Material	Nickel Plated OFHC Copper
Attachment	Flanged
Optical Access	0, 2, or 4 (customer specified)

Cryostat Footprint

Overall Length	603 mm (23.72 in)
Motor Housing Diameter	114 mm (4.5 in)
Rotational Clearance	200 mm (8 in) with "G" Configuration


Cryocooler Model

		DE-202AI		DE-202A(T)I		DE-202PI		DE-202SI	
	Frequency	60 Hz	50 Hz	60 Hz	50 Hz	60 Hz	50 Hz	60 Hz	50 Hz
Base Temperature		<9K	<9K	<9K	<9K	<5.5K	<5.5K	<4.2K	<4.2K
Cooling Capacity*	4.2K	-	-	-	-	-	-	0.1W	0.08W
	10K	0.5W	0.4W	0.7W	0.56W	1W	0.8W	1.2W	1W
	20K	2.5W	2W	3.7W	3W	3.5W	2.8W	4W	3.2W
	77K	4W	3.2W	6W	4.8W	3.5W	2.8W	4W	3.2W
Radiation Shield Cooling Capacity		10W	8W	15W	12W	10W	8W	10W	8W
Cooldown Time	20K	50 min	60 min	35 min	42 min	60 min	72 min	60 min	72 min
	Base Temperature	70 min	84 min	50 min	60 min	90 min	108 min	90 min	108 min
Compressor Model		ARS-4HW		ARS-4HW		ARS-4HW		ARS-4HW	
Typical Maintenance Cycle		12,000 hours		12,000 hours		12,000 hours		12,000 hours	



Optical Cryostat - Large Sample Space


CS202*I-DMX-4SS Outline Drawing

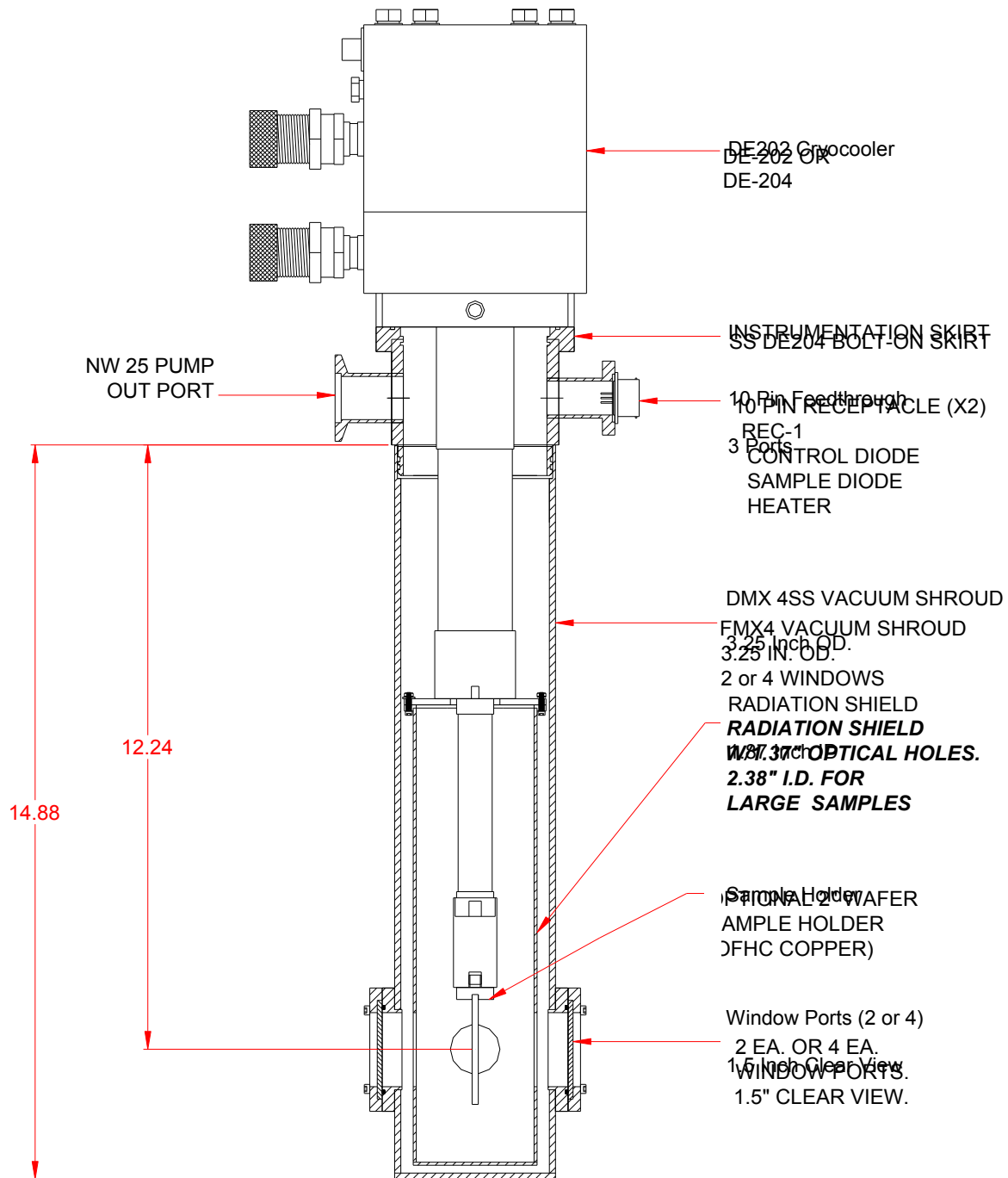




Optical Cryostat - Large Sample Space

Sample Space

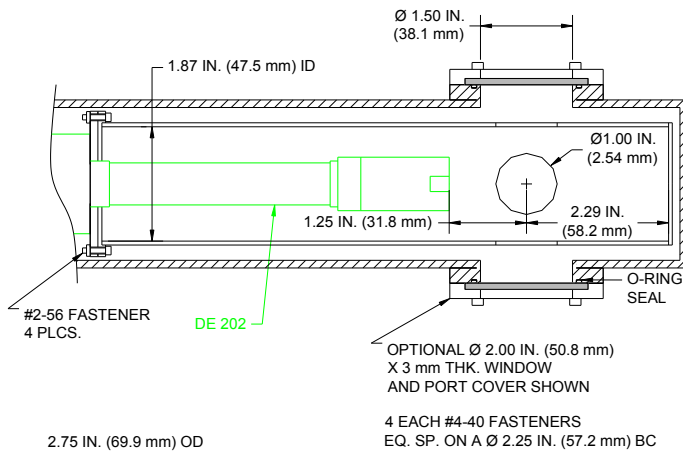

ARS-4HW Compressor

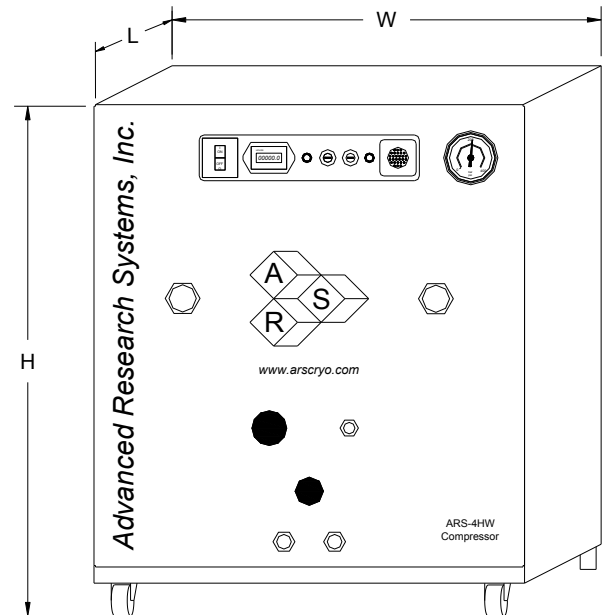

Ø2.50 IN. (63.5 mm)

Compressor Model

ARS-4HW

12.25

14.88

		Frequency		60 Hz	50 Hz
Standard Voltage	Min			208 V	190 V
	Max			230 V	210 V
Transformer Options	10%			220 V, 230 V	
	15%			240 V	
Power Usage	Single Phase			3.6 kW	3.0 kW
Refrigerant Gas				99.999% Helium Gas, Pre-Charged	
Noise Level				60 dBA	
Ambient Temperature				12 - 40 C (54 - 104 F)	
Cooling Water	Consumption			2.3 L / min (0.6 Gal. / min)	
	Temperature			10 - 35 C (50-95 F)	
	Connection			3/8 in. Swagelok Fitting	
Dimensions:	L			483 mm (19 in)	
	W			434 mm (17.1 in)	
	H			516 mm (20.3 in)	
Weight				72 kg (160 lbs)	
Typical Maintenance Cycle				12,000 hours	
Water Recirculation Option				CoolPac Compatible	